

Clinet® Anti-Algae

Liquid for inhibiting algae growth



Efficiency for lab baths

- For water baths hygiene
- Reduces the water consumption
- Change the water every month only
- Easy-to-use thanks to its designed bottle

Clinet® Anti-Algae maintains and cleans the water baths and ultrasonic trays.

interscience

Applications

Prevents algae for water baths, pool baths...

How to use

- Use it diluted at 0.1 %:
1 dose of 20 mL for 20 L of demineralised water
- Renew the necessary dose every week
- Change the water bath every month

Composition

Ammonium quaternary, calcium, preservative

Precautions

- Use exclusively demineralized or deionized water because the algicide action is diminished by hard water.
- Do not mix with anion-active detergents and/or strong oxydants (such as bleach...)
- In case of contact with the eyes, rinse abundantly with water and consult a specialist
- Do not swallow
- Keep out of reach of children



For more info, the safety data sheet is available on request.

Technical specifications

Clinet® Anti-Algae	
Reference	310 400
Box of	1 L
Size	34 x 25.5 x 18.5 cm
Gross weight	1.1 kg

Clinet® range



Clinet® 101
Disinfecting wipes
Ref.: 310 101




Clinet® 202
Disinfecting wipes
in individual pack
Ref.: 310 202



Clinet® 303
Ready-to-use
disinfecting spray
Ref.: 310 303

Discover our complete range for microbiology

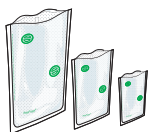

BagFilter®
Sterile filter bags



DiluFlow®
Gravimetric dilutors



BagMixer®
Lab blenders



FlexiPump®
Peristaltic pumps



easySpiral Dilute®
Automatic diluter and platers



Scan®
Automatic colony counters
Inhibition zone readers

Distributor

interscience INTERNATIONAL

30, chemin du Bois des Arpents - 78860 St Nom - FRANCE
T: +33 (0)1 34 62 62 61 - F: +33 (0)1 34 62 43 03 - E-mail: info@interscience.com

interscience USA & CANADA

199 Weymouth street - ROCKLAND - MA 02370 - USA
T: +1 781 792 2133 - F: +1 781 792 2134 - E-mail: sales.usa@interscience.com

interscience CHINA

上海市徐汇区吴兴路277号锦都大厦718室 - 200030
电话: +86 (0)21-64739390 - 手机: +86 189 3097 0733 - 邮址: sales.china@interscience.com

interscience SOUTH-EAST ASIA

541 Orchard road - 09-01 Liat Towers - SINGAPORE 238881
T: +65 6933 1389 - F: +65 8549 1217 - E-mail: sales.asia@interscience.com

www.interscience.com