

Thermo Scientific Finnpipette F1 Pipette

Thermo Scientific™ Finnpipette™ F1 pipettes deliver performance with outstanding precision in a lightweight design that maximizes user comfort and reduces strain. It's the ideal system for applications using expensive reagents and valuable samples, thanks to its super blow-out feature for effective delivery of microvolumes, along with a secure volume-lock capability that prevents unintended volume shift during operation.

Ergonomically In-touch

Lightweight handle design combined with low plunger, tip attachment and ejection forces help reduce the risk of Repetitive Strain Injury (RSI).

Silent, Effortless Volume Adjustment

Enjoy the new silent and lighter volume adjustment mechanism with over 50% lower forces compared to earlier models. Improved grip on the volume adjustment knob ensures effortless and precise adjustment.

Secure Volume Lock

The secure volume setting prevents undesirable volume drift during operation. Simply lift the plunger up to set volume and press down to lock in place with minimum effort.

Adjustable Comfort

The 120° adjustable finger rest – for both right- and left-handed operator comfort – allows the hand to relax during pipetting cycles.

Super Blow-out

150% increase in air boost to ensure efficient delivery of micro-volumes and prevent capillary action in 50 µL models and below.

* Validity may vary by country.



Wide Selection of Finntip Pipette Tips

Finntips are designed and manufactured expressly for Finnpipette models, enabling optimal performance, precision and accuracy.

5-Year Warranty

With convenient online registration*.



The 120° adjustable finger rest

Thermo Scientific Finnpiptette F1 Pipettes

Item No.	Color Key/Description	Increment	Volume (μL)	Inaccuracy		Imprecision		Compatible Tip
				μL	%	s.d.* μL	cv%	
Finnpiptette F1 Single Channel Pipettes								
4641010N	● Finnpiptette F1, 0.2-2 μL , micro	0.002 μL	2	± 0.050	± 2.50	0.04	2.00	Finntip 10, Flex 10
			1	± 0.040	± 4.00	0.035	3.50	
			0.2	± 0.024	± 12.00	0.02	10.00	
4641020N	● Finnpiptette F1, 0.5-5 μL , micro	0.01 μL	5	± 0.075	± 1.50	0.05	1.00	Finntip 10, Flex 10
			2.5	± 0.063	± 2.50	0.038	1.50	
			0.5	± 0.030	± 6.00	0.025	5.00	
4641030N	● Finnpiptette F1, 1-10 μL , micro	0.02 μL	10	± 0.100	± 1.00	0.05	0.50	Finntip 10, Flex 10
			5	± 0.075	± 1.50	0.04	0.80	
			1	± 0.025	± 2.50	0.02	2.00	
4641040N	● Finnpiptette F1, 1-10 μL	0.02 μL	10	± 0.100	± 1.00	0.08	0.80	Finntip 250, Flex 200
			5	± 0.075	± 1.50	0.04	0.80	
			1	± 0.035	± 3.50	0.03	3.00	
4641050N	● Finnpiptette F1, 2-20 μL , micro	0.02 μL	20	± 0.20	± 1.00	0.08	0.40	Finntip 50
			10	± 0.15	± 1.50	0.06	0.60	
			2	± 0.06	± 3.00	0.05	2.50	
4641060N	● Finnpiptette F1, 2-20 μL	0.02 μL	20	± 0.20	± 1.00	0.08	0.40	Finntip 250, Flex 200
			10	± 0.15	± 1.50	0.06	0.60	
			2	± 0.06	± 3.00	0.05	2.50	
4641130N	● Finnpiptette F1, 5-50 μL , micro	0.1 μL	50	± 0.30	± 0.60	0.15	0.30	Finntip 50
			25	± 0.25	± 1.00	0.13	0.50	
			5	± 0.15	± 3.00	0.125	2.50	
4641140N	● Finnpiptette F1, 5-50 μL	0.1 μL	50	± 0.30	± 0.60	0.15	0.30	Finntip 250, Flex 200
			25	± 0.25	± 1.00	0.13	0.50	
			5	± 0.15	± 3.00	0.125	2.50	
4641070N	● Finnpiptette F1, 10-100 μL	0.2 μL	100	± 0.80	± 0.80	0.20	0.20	Finntip 250, Flex 200
			50	± 0.60	± 1.20	0.20	0.40	
			10	± 0.30	± 3.00	0.10	1.00	
4641080N	● Finnpiptette F1, 20-200 μL	0.2 μL	200	± 1.2	± 0.60	0.40	0.20	Finntip 250, Flex 200
			100	± 1.0	± 1.00	0.40	0.40	
			20	± 0.36	± 1.80	0.14	0.70	
4641090N	● Finnpiptette F1, 30-300 μL	1 μL	300	± 1.8	± 0.60	0.60	0.20	Finntip 300, Flex 300
			150	± 1.5	± 1.00	0.60	0.40	
			30	± 0.45	± 1.50	0.18	0.60	
4641100N	● Finnpiptette F1, 100-1000 μL	1 μL	1000	± 6.0	± 0.60	2.00	0.20	Finntip 1000, Flex 1000
			500	± 4.0	± 0.80	1.50	0.30	
			100	± 1.0	± 1.00	0.60	0.60	
4641110N	● Finnpiptette F1, 0.5-5 mL	0.01 mL	5000	± 25.0	± 0.50	10.00	0.20	Finntip 5 mL
			2500	± 17.5	± 0.70	7.50	0.30	
			500	± 10.0	± 2.00	4.00	0.80	
4641120	● Finnpiptette F1, 1-10 mL	0.02 mL	10000	± 50.0	± 0.50	20.00	0.20	Finntip 10 mL, Flex 10ml Ext
			5000	± 40.0	± 0.80	15.00	0.30	
			1000	± 20.0	± 2.00	8.00	0.80	
Finnpiptette F1 Multichannel Pipettes								
4661000N	● Finnpiptette F1, 8-ch, 1-10 μL	0.02 μL	10	± 0.240	± 2.40	0.16	1.60	Finntip 10, Flex 10
			5	± 0.200	± 4.00	0.15	3.00	
			1	± 0.120	± 12.00	0.08	8.00	
4661010N	● Finnpiptette F1, 8-ch, 5-50 μL	0.1 μL	50	± 0.75	± 1.50	0.35	0.70	Finntip 250, Flex 200
			25	± 0.625	± 2.50	0.30	1.20	
			5	± 0.25	± 5.00	0.10	2.00	
4661020N	● Finnpiptette F1, 8-ch, 10-100 μL	0.2 μL	100	± 1.30	± 1.30	0.50	0.50	Finntip 250, Flex 200
			50	± 1.25	± 2.50	0.60	1.20	
			10	± 0.50	± 5.00	0.20	2.00	
4661030N	● Finnpiptette F1, 8-ch, 30-300 μL	1 μL	300	± 3.0	± 1.00	0.90	0.30	Finntip 300, Flex 300
			150	± 2.25	± 1.50	0.75	0.50	
			30	± 1.5	± 5.00	0.60	2.00	
4661040N	● Finnpiptette F1, 12-ch, 1-10 μL	0.02 μL	10	± 0.240	± 2.40	0.16	1.60	Finntip 10, Flex 10
			5	± 0.200	± 4.00	0.15	3.00	
			1	± 0.120	± 12.00	0.08	8.00	
4661050N	● Finnpiptette F1, 12-ch, 5-50 μL	0.1 μL	50	± 0.75	± 1.50	0.35	0.70	Finntip 250, Flex 200
			25	± 0.625	± 2.50	0.30	1.20	
			5	± 0.25	± 5.00	0.10	2.00	
4661060N	● Finnpiptette F1, 12-ch, 10-100 μL	0.2 μL	100	± 1.30	± 1.30	0.50	0.50	Finntip 250, Flex 200
			50	± 1.25	± 2.50	0.60	1.20	
			10	± 0.50	± 5.00	0.20	2.00	
4661070N	● Finnpiptette F1, 12-ch, 30-300 μL	1 μL	300	± 3.0	± 1.00	0.90	0.30	Finntip 300, Flex 300
			150	± 2.25	± 1.50	0.75	0.50	
			30	± 1.5	± 5.00	0.60	2.00	
4661080N	● Finnpiptette F1, 16-ch, 1-10 μL	0.02 μL	10	± 0.240	± 2.40	0.16	1.60	Finntip 20
			5	± 0.200	± 4.00	0.15	3.00	
			1	± 0.120	± 12.00	0.08	8.00	
4661090N	● Finnpiptette F1, 16-ch, 5-50 μL	0.1 μL	50	± 0.75	± 1.50	0.35	0.70	Finntip 50
			25	± 0.625	± 2.50	0.30	1.20	
			5	± 0.25	± 5.00	0.10	2.00	

*Standard Deviation

Finnpipette F1 Fixed Volume Single Channel Pipettes

Item No.	Description	Volume (μ L)	Inaccuracy		Imprecision		Compatible Tip
			μ L	%	s.d.* μ L	cv%	
4651020N	Finnpipette F1 Fixed 10 μ L	10	± 0.090	± 0.90	0.08	0.80	Finntip 250, Flex 200
4651130N	Finnpipette F1 Fixed 20 μ L	20	± 0.14	± 0.70	0.10	0.50	Finntip 250, Flex 200
4651030N	Finnpipette F1 Fixed 25 μ L	25	± 0.15	± 0.60	0.125	0.50	Finntip 250, Flex 200
4651040N	Finnpipette F1 Fixed 50 μ L	50	± 0.30	± 0.60	0.20	0.40	Finntip 250, Flex 200
4651050N	Finnpipette F1 Fixed 100 μ L	100	± 0.40	± 0.40	0.30	0.30	Finntip 250, Flex 200
4651140N	Finnpipette F1 Fixed 200 μ L	200	± 0.80	± 0.40	0.60	0.30	Finntip 250, Flex 200
4651060N	Finnpipette F1 Fixed 250 μ L	250	± 1.0	± 0.40	0.75	0.30	Finntip 1000, Flex 1000
4651070N	Finnpipette F1 Fixed 500 μ L	500	± 1.5	± 0.30	1.50	0.30	Finntip 1000, Flex 1000
4651080N	Finnpipette F1 Fixed 1000 μ L	1000	± 3.0	± 0.30	3.00	0.30	Finntip 1000, Flex 1000
4651090N	Finnpipette F1 Fixed 2000 μ L	2000	± 6.0	± 0.30	4.00	0.20	Finntip 5 mL
4651100N	Finnpipette F1 Fixed 3000 μ L	3000	± 9.0	± 0.30	6.00	0.20	Finntip 5 mL
4651110N	Finnpipette F1 Fixed 5000 μ L	5000	± 15.0	± 0.30	10.0	0.20	Finntip 5 mL
4651120N	Finnpipette F1 Fixed,10000 μ L	10000	± 30.0	± 0.30	20.0	0.20	Finntip 10 mL, Flex 10 mL Ext

*Standard Deviation

Finnpipette F1 Good Laboratory Pipetting Kits

Item No.	Description	Kit Includes
4700850N	Finnpipette F1 GLP Kit 1	Finnpipette F1 1-10 μ L single channel Finnpipette F1 10-100 μ L single channel Finnpipette F1 100-1000 μ L single channel F-Series Stand 3 Racks of Finntip Flex tips
4700860N	Finnpipette F1 GLP Kit 2	Finnpipette F1 0.2-2 μ L single channel Finnpipette F1 2-20 and 20-200 μ L single channels Finnpipette F1 100-1000 μ L single channel F-Series Stand 4 Racks of Finntip/Finntip Flex tips
4700865N	Finnpipette F1 GLP Kit 3	Finnpipette F1 10-100 μ L single channel Finnpipette F1 100-1000 μ L single channel Finnpipette F1 1-10 mL single channel F-Series Stand 3 Racks of Finntip/Finntip Flex tips
4701060N	Finnpipette F1 GLP Kit 4	Finnpipette F1 2-20 μ L single channel Finnpipette F1 20-200 μ L single channel Finnpipette F1 100-1000 μ L single channel F-Series Stand 3 Racks of Finntip Flex tips



Remember to Register Your Pipette
for a **Free** Extended Warranty!
www.thermoscientific.com/pipetwarranty

Learn more at thermoscientific.com/finnpipette

© 2016 Thermo Fisher Scientific Inc. All rights reserved.

All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Please consult your local sales representative for details.

Australia +61 39757 4300

Austria +43 1 801 40 0

Belgium +32 53 73 42 41

China +86 21 6865 4588 or

+86 10 8419 3588

France +33 2 2803 2180

Germany national Toll-Free 0800 1 536 376

Germany International +49 6184 90 6000

India Toll-Free 1800 22 8374

India +91 22 6716 2200

Italy +39 02 95059 554

Japan +81 45 453 9220

Netherlands +31 76 579 55 55

New Zealand +64 9 980 6700

Nordic/Baltic/CIS Countries

+358 9 329 10200

Russia +7 812 703 42 15

Spain/Portugal +34 93 223 09 18

Switzerland +41 44 454 12 22

UK/Ireland +44 870 609 9203

USA/Canada +1 866 984 3766

Other Asian countries +852 2885 4613

Countries Not Listed +49 6184 90 6000

SSHPFinnF10716
1517680

ThermoFisher
SCIENTIFIC